

Tryptic Standard I Nutrient Agar 90 mm CS

Technical Data Sheet													
Use in	<ul style="list-style-type: none"> • Pharmaceutical industry in clean rooms and isolators • For industrial, laboratory & research applications only • Basic medium according to Ph. Eur. and USP 												
Use for	<ul style="list-style-type: none"> • Detection of aerobic and anaerobic micro-organisms • Active and passive air sampling, personnel monitoring • Isolation and growth of fastidious bacteria, yeasts and moulds • Neutralization of residues of disinfectants 												
Typical composition per liter	<table style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 30%;">Peptone</td> <td style="width: 10%;">15 g</td> <td style="width: 30%;">Glucose x H₂O</td> <td style="width: 30%;">1.1 g</td> </tr> <tr> <td>Yeast extract</td> <td>3 g</td> <td>Agar</td> <td>15 g</td> </tr> <tr> <td>NaCl</td> <td>6 g</td> <td></td> <td></td> </tr> </table> <p>This medium can be adjusted / or supplemented according to the performance criteria required.</p>	Peptone	15 g	Glucose x H ₂ O	1.1 g	Yeast extract	3 g	Agar	15 g	NaCl	6 g		
Peptone	15 g	Glucose x H ₂ O	1.1 g										
Yeast extract	3 g	Agar	15 g										
NaCl	6 g												
Irradiation	<ul style="list-style-type: none"> • Not irradiated 												
Filling volume	<ul style="list-style-type: none"> • 23-26 mL 												
Packaging	<ul style="list-style-type: none"> • Single bagged, staples of 10 plates • Transparent • High barrier foil against desiccation • 6 staples of 10 plates per packaging unit • Temperature insulated handle-bag in the cardboard-boxes 												
Plates per box	<ul style="list-style-type: none"> • 60 plates (6 staples with 10 plates each) 												
Shelf life	<ul style="list-style-type: none"> • 8 months from production date – preliminary expiry date 												
Storage conditions	<ul style="list-style-type: none"> • Recommended storage temperature: 15-25 °C • Should be stored at temperatures as stable as possible • Store protected from light exposure • Before use: it is recommended to keep the plates upright (agar on the lower part, lid on the upper part) to avoid formation of extra condensation • After use: it is recommended to keep the plates upside down (agar on the upper part, lid on the lower part) to reduce the risk of accumulation of condensation during incubation which can affect colony formation 												
Label	<ul style="list-style-type: none"> • On the side of the bottom part of the dish 												
Label information	<ul style="list-style-type: none"> • Product name: Std.-I • Expiry date: YYYYMMDD → MMM in letters (e.g.: 2026Nov04) • Lot-number • Individual number • Barcode 												

Technical Data Sheet	
Barcode	<ul style="list-style-type: none"> • 2-dimensional (data matrix), 20 digits: • Digits 1-3: Art.-No. • Digits 4-9: Lot-Number • Digits 10-14: Individual-Number • Digits 15-20: Date (YYMMDD)
Delivery	<ul style="list-style-type: none"> • Temperature controlled delivery on request • For shipments of larger amounts plastic pallets in Euro-size can be used
Petri dish	<ul style="list-style-type: none"> • Locking lid 90 mm plate, made from polystyrene • Long incubations possible – due to high filling volume • Long expositions possible – due to specific design of plate • Incubations in vent and closed position possible
Lid positions	<ul style="list-style-type: none"> • All plates are delivered in the non-locked position • The plate contains two locked positions. If turning the lid clockwise the locked positions are in the following order: <ol style="list-style-type: none"> 1. Vent position 2. Closed position
Aerobic incubation	<ul style="list-style-type: none"> • The closed position provides ideal incubation conditions for aerobic microorganisms and limits the dehydration of the agar during incubation • For long incubation of aerobic microorganisms, the closed position is recommended • To lock the lid in the closed position, turn the lid clockwise into the final stop position
Anaerobic incubation	<ul style="list-style-type: none"> • The vent position is ideal for anaerobic incubations, as it allows an easy and effective removal of oxygen under anaerobic incubation conditions • Incubate in anaerobic incubator, anaerobic jar or suitable equipment <ol style="list-style-type: none"> 1. First option: <ul style="list-style-type: none"> • Turn the lid clockwise into the final stop position • Turn the lid one click counterclockwise to the vent position 2. Second option: <ul style="list-style-type: none"> • Turn the lid clockwise directly into the first locked position
Place of production	PharmaMedia Dr. Müller GmbH Gustav-Throm-Str. 1, 69181 Leimen - Germany

Quality control, Certificates																					
Certificates	Each lot of product can be obtained with a certificate of analysis (CoA):																				
	<table border="1"> <thead> <tr> <th colspan="2">Physico-chemical test parameters:</th> </tr> </thead> <tbody> <tr> <td>Appearance</td> <td>Clear, yellowish</td> </tr> <tr> <td>pH value</td> <td>7.3 – 7.7</td> </tr> <tr> <td>Filling volume</td> <td>23 – 26 mL</td> </tr> </tbody> </table>	Physico-chemical test parameters:		Appearance	Clear, yellowish	pH value	7.3 – 7.7	Filling volume	23 – 26 mL												
	Physico-chemical test parameters:																				
	Appearance	Clear, yellowish																			
	pH value	7.3 – 7.7																			
	Filling volume	23 – 26 mL																			
	<table border="1"> <thead> <tr> <th colspan="5">Growth Promotion test: 10-100 CFU</th> </tr> </thead> <tbody> <tr> <td><i>E. coli</i></td> <td>ATCC 8739</td> <td>36-38 °C</td> <td>1 day</td> <td>50-200%</td> </tr> <tr> <td><i>S. Thyphimurium</i></td> <td>ATCC 14028</td> <td>36-38 °C</td> <td>1 day</td> <td>50-200%</td> </tr> <tr> <td><i>S. Abony</i></td> <td>NCTC 6017</td> <td>36-38 °C</td> <td>1 day</td> <td>50-200%</td> </tr> </tbody> </table>	Growth Promotion test: 10-100 CFU					<i>E. coli</i>	ATCC 8739	36-38 °C	1 day	50-200%	<i>S. Thyphimurium</i>	ATCC 14028	36-38 °C	1 day	50-200%	<i>S. Abony</i>	NCTC 6017	36-38 °C	1 day	50-200%
	Growth Promotion test: 10-100 CFU																				
	<i>E. coli</i>	ATCC 8739	36-38 °C	1 day	50-200%																
	<i>S. Thyphimurium</i>	ATCC 14028	36-38 °C	1 day	50-200%																
<i>S. Abony</i>	NCTC 6017	36-38 °C	1 day	50-200%																	
<table border="1"> <thead> <tr> <th colspan="2">Sterility control</th> </tr> </thead> <tbody> <tr> <td></td> <td>Conform</td> </tr> </tbody> </table>	Sterility control			Conform																	
Sterility control																					
	Conform																				
Certificate of origin	All media lots produced by PMM can be obtained with a Certificate of Origin (CoO). All animal derived raw materials are specified as follows:																				
	<ul style="list-style-type: none"> • Raw material • Tissue • Animal source • Country of origin • Infectivity category (acc. to TSE guideline: EMA/410/01 current version) 																				
BSE policy	<ul style="list-style-type: none"> • In compliance with the current note for guidance on minimizing the risk of transmitting animal spongiform encephalopathy via human or veterinary medicinal products, we check the CoO of raw material in respect to the specified animal source, the country of origin and the infectivity category. We neither store or process ruminant raw materials obtained from high infectivity tissues (IA) nor ruminant raw materials whose animal source originates from countries or regions with an undetermined risk (cat C/GBR IV). 																				

Safety Data	
Toxic ingredients	<ul style="list-style-type: none"> • None
Basic composition	<ul style="list-style-type: none"> • See typical composition
Solvent content	<ul style="list-style-type: none"> • None
Safety data sheet required	<ul style="list-style-type: none"> • Not mandatorily required